




Unpacking Instructions

GX-281 Liquid Handler

The GX-281 Liquid Handler is delivered with most major components already assembled. Keep the original container and packing assembly in case the GX-281 Liquid Handler must be returned to the factory.

 At least two people should lift the liquid handler off the base of the packing container, as it weighs over 68 kg (150 lbs).

- 1 Remove the box containing the shield (if included) from the top of the liquid handler.
- 2 Remove the accessory boxes from the locator plate of the liquid handler.
- 3 Lift the liquid handler off its base platform and place it on a lab bench or cart.

To lift the liquid handler:

- a) Grip the liquid handler under the base plate.
- b) Lift the unit up and out of the foam packing material.

NOTICE: Do not attempt to lift the instrument from the Y-arm (the horizontal arm). Always lift the instrument from its base.

After the unit has been unpacked, you should have the following items:

- GX-281 Liquid Handler
- Z-arm
- GX-281 accessory package
- GX Prep Solvent System (included with the GX-281 Prep Liquid Handler)
- GX Analytical Solvent System (included with the GX-281 Analytical Liquid Handler)
- Power cord, 220V
- Power cord, 110V
- GX-281 Offset Utility Kit
- Doc CD, GX-281 Liquid Handler
- Items Included Checklist, GX-281
- Declaration of Conformity GX-281
- Unpacking the GX-281 Liquid Handler
- IQ Procedure GX-281 Liquid Handler
- QC Checklist GX-281 Liquid Handler

