

# INSTALLING YOUR SINGLE™ PIPETTE HOLDER

---

*for PIPETMAN®, MICROMAN® and  
DISTRIMAN®*

Quick and easy **adhesive attachment** for use on any smooth clean, dry surface or **screw fitting** for use on rough surfaces.

*There are two ways of fixing your SINGLE to a wall, according to the nature of the surface.*

## **For fixing on a smooth surface – use the double-sided adhesive tape.**

The surface must be clean, degreased, and dry. The ideal temperature for applying the tape is between 21°C and 38°C.

- 1) First, clean the surface with an ethanol and water based solvent. You are advised to follow the safety instructions given by the supplier when you use the solvent.
- 2) Remove the backing from one side of the tape and carefully stick the tape onto the back of your SINGLE.

- 3) Remove the backing from the other side of the tape, place the SINGLE at the required position and press it hard against the surface, pushing through the inside of the SINGLE.
- 4) Wait 24 hours before using the SINGLE.

**For fixing on a rough surface - use the wall-plugs and screws**

- 1) Drill two holes at the appropriate position using a 6-mm drill.
- 2) Push a wall-plug into each of the two holes.
- 3) Screw in the two screws, leaving a 2-mm gap between the heads of the screws and the wall.
- 4) Position slots in the back of the SINGLE over the screws and push down to install it. Tighten the screws, if necessary, to maintain the SINGLE firmly against the wall.

ORDERING REF. N° F161406

LT801302/D - Dec. 2016